

# Technical Data Sheet

## CellScrew® mini TriDock



*Document version number:* ..... 2.0

*Date of issue:* ..... July 24, 2024



## Index

- 1. Product description . . . . . 3
- 2. Technical overview . . . . . 3
- 3. Instructions for use . . . . . 4
- 4. Technical support . . . . . 5



## 1. Product description

The CellScrew® mini TriDock is a reusable device designed for use with the CellScrew® mini, enabling the simultaneous operation of three CellScrew® minis. It is manufactured in an additive manufacturing process and is made of plant-based polylactic acid (PLA). The CellScrew® mini TriDock has the same outer diameter as the CellScrew® 6K and 10K, making it compatible with standard roller devices. Key data is provided below.

## 2. Technical overview

### Dimensions

Dimension	Unit	Value
Diameter	[mm]	121
Total height	[mm]	192
Bottom part height	[mm]	169
Lid height	[mm]	60
Screw height	[mm]	32


### Properties

<b>Description:</b>	Green Elephant Biotech GmbH CellScrew® mini TriDock GEB-CS-mini-TriDock
<b>Packaging dimensions:</b>	130 mm x 130 mm x 300 mm
<b>Weight:</b>	190 g
<b>Material:</b>	Polylactic acid
<b>Color:</b>	White
<b>Temperature range:</b>	- 80 °C to 45 °C
<b>Storage:</b>	Room temperature, sunlight protected
<b>Centrifugability:</b>	No
<b>Autoclavability:</b>	No
<b>Recommended rotational speed:</b>	0.2 – 2 rpm
<b>Packing unit:</b>	1 CellScrew® mini TriDock, 1 piece per box
<b>Quality management system:</b>	DIN EN ISO 9001:2015



## 3. Instructions for use

The CellScrew® mini TriDock consists of three components: a bottom part, a lid, and a screw. The screw secures the lid to the central bar of the bottom part.

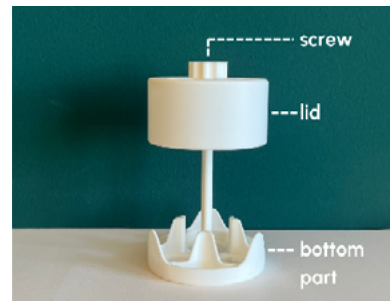

Three CellScrew® minis fit into the CellScrew® mini TriDock. Place them in the designated cut-outs of the bottom part.



Align the round recesses in the lid with the caps and filters of the CellScrew® minis.



Close the lid by tightening the screw.



The CellScrew® mini TriDock is operated similarly to a CellScrew® 6K or 10K, by placing it on an angled (10° - 15°) roller device. The CellScrew® mini TriDock must rotate clockwise when observed from above, with a rotational speed ranging from 0.2 to 2 rpm.





Remove the CellScrew® mini from the CellScrew® mini TriDock for cell inoculation and harvesting steps, when working inside an asptic working bench.

Inoculate the CellScrew® mini with the desired cell concentration and start incubation using the CellScrew® mini TriDock. Put it on an angled (10 ° - 15 °) roller device and rotate it clockwise at 0.2 - 2 rpm.

After cell expansion remove the culture medium and add your washing solution. Wash the cells by holding the CellScrew® mini at an angle and rotating it clockwise by hand. Remove the washing solution and add trypsin or another detachment agent to the CellScrew® mini. Distribute the detachment agent by holding the CellScrew® mini at an angle and rotating it clockwise by hand. Incubate for the desired detachment time at the desired temperature using the CellScrew® mini TriDock. Put it on an angled (10 ° - 15 °) roller device and rotate it clockwise at 0.2 - 2 rpm during detachment.

Stop the detachment reaction with FBS containing medium or a different suitable stopping reagent. Hold the CellScrew® mini at an angle and tap it firmly, while rotating it clockwise by hand to detach the cells. Bring the CellScrew® mini to an upright position and transfer the cell suspension into a harvest vessel for further use.

Detailed instructions are available in the [Technical data sheet](#) for the CellScrew® mini.

## 4. Technical support

For technical support, questions or remarks please contact the technical department of Green Elephant Biotech GmbH via e-mail ([tech-service@greenebt.com](mailto:tech-service@greenebt.com)).